



ND-1000 CALIBRATION CHECK

INSTALL SOFTWARE

Download and install the latest version of the ND-1000 Calibration Check Software from the Support Section on the NanoDrop web site at www.nanodrop.com.

CALIBRATION PROCEDURE

1. Open the ND-1000 Calibration Check Software and follow the prompts in the Customer Guidance section of the software.
2. Before opening the ampule of CF-1 Calibration Fluid, shake vigorously to ensure solution is thoroughly mixed.
3. Tap upright ampule lightly to ensure all solution is collected in the bottom portion of the ampule.
4. Carefully break neck of ampule to open CF-1 Calibration Fluid.
5. If average value is unacceptable, contact NanoDrop Technologies @ 302-479-7707.

NOTE: The CF-1 Calibration fluid is intended for single use. Exposure to the environment or transferring the fluid to another container can cause a significant concentration change.

